**Chromatography** Chromatographie

**Multidimensional  
Chromatography** Multidimensionale  
Chromatographie

# Multidimensional Chromatography

Multidimensional methods allow the analysis of complex samples in a few steps. Beside the coupling of several columns with the same properties for highest resolution, combinations of chemically different stationary phases are also used. These approaches enable the increase of the peak capacity and can help to avoid co-elution as can happen in a single dimension chromatography.

In HPLC and GC practice different concepts are used and common:

- peak cutting and the selective transfer of peaks or chromatogram sections onto a second column (heart-cut method)
- continuous two-dimensional chromatography (comprehensive chromatography).

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#### Comparison between Traditional Deans Switching and Multi-Deans Switching

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### Mechanism of Multi-Deans Switching

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$P - \Delta P2 > P - \Delta P1$  1. DET

# Heart-Cut Gas Chromatography/Gas-Chromatographie

The multidimensional gas chromatography adds a third analytical dimension. With the new Multi-Deans Switch technology the operator can define as many heart-cuts to be transferred to the second column as needed for his target compound analysis.

The MDGC-2010 system can be successfully employed for a wide range of applications in, e.g. flavors and fragrances, environmental, food and clinical analysis. It sets a new standard in multidimensional gas chromatography. The Multi-Deans Switching offers all possibilities of the heart-cut technique without its previous drawbacks.

- very easy method development: just three steps are needed to set up a multidimensional method
- as many heart-cuts as required without any shifts of the very precise retention times
- very easy quantitation
- all operation parameters are controlled via MDGCsolution software
- for detection all conventional GC detectors as well as the extremely sensitive and fast Quadrupole GCMS-QP2010 Plus can be used

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## LC x LC

Comprehensive LC (also called LC x LC) couples two liquid chromatography based separation methods and enables the separation of complex samples as well as the identification of co-eluting peaks. Different separation mechanisms should be chosen in order to achieve maximum possible resolution.

- two binary gradient systems enable maximum flexibility in choosing the separation method and mechanism
- detection after the second dimension only minimizes the transfer volume between the separation steps
- for the sample transfer a switching valve with two sample loops is used and operated in continuous mode
- all types of LC detectors and mass spectrometers can be used

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## GC x GC

### GC x GC (qMS)

Comprehensive chromatography is a true breakthrough in qualitative and quantitative high-end analysis for samples of natural origin from e.g. petrochemical, cosmetic, flavors and fragrances and food industry. It is based on two columns of different polarities.

The key component of GC x GC (qMS) is the patented Zoex modulator connecting both columns. It focuses the compounds from the first column and injects them into the second column.

Correct representation of the very narrow peaks obtained by:

- fast GC-detectors with a minimum filter time constant of 4 ms and a maximum data acquisition frequency of 250 Hz (see chromatography-brochure GC/GCMS)
- GCMS-QP2010 Plus featuring the fastest acquisition rates with a maximum of 50 scans per second

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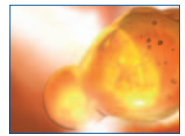
## LC-GC coupling

For complex samples it may not be sufficient to couple two GC columns of different polarity to characterize all components. In this case, LC can be coupled to GC. Whereas LC separates compounds by their polarity or size, compounds in GC using non-polar columns are separated by their boiling points.

These different separation mechanisms can be used for the separation of co-elutions in various applications, e.g. characterization of olive oil or polymers or the determination of pesticides in complex matrices. Transfer of the fractions to the GC is performed via a special syringe using the AOC-5000 autosampler.

In principle, LC and GC can be coupled in two different ways:

- inject LC fractions in pre-set time intervals into the GC system (comprehensive chromatography)
- selected parts of the LC chromatogram are injected into the GC for further analysis (heart-cut method).

In the latter case LC serves as sample preparation for the GC. For detection all conventional GC detectors can be used. For correct identification and quantitation of the compounds in the usually very complex samples, however, in most cases a Quadrupole GCMS system is used.

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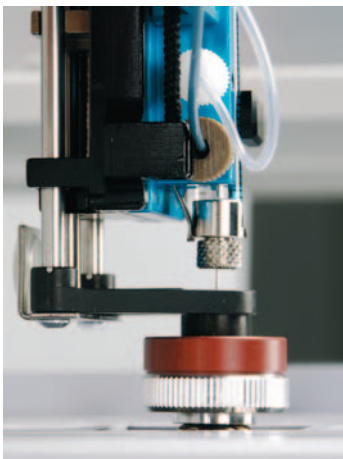
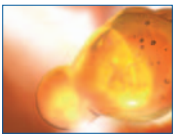
## Data acquisition and evaluation software packages

### LabSolutions software

- acquires and processes data in multidimensional and comprehensive chromatography

### ChromSquare software

- easy creation of contour plots for further data evaluation in comprehensive chromatography
- identification of compounds in GC-MS applications directly from the contour plot using library searching with linear retention index search

### GCImage software (Zoex)

- for comprehensive GC (GC x GC [qMS]) this well-established tool is perfect for qualitative and quantitative data analysis

### Unique LC x GC software

- controls the complete LC and GC systems including the AOC-5000 autosampler from one user interface

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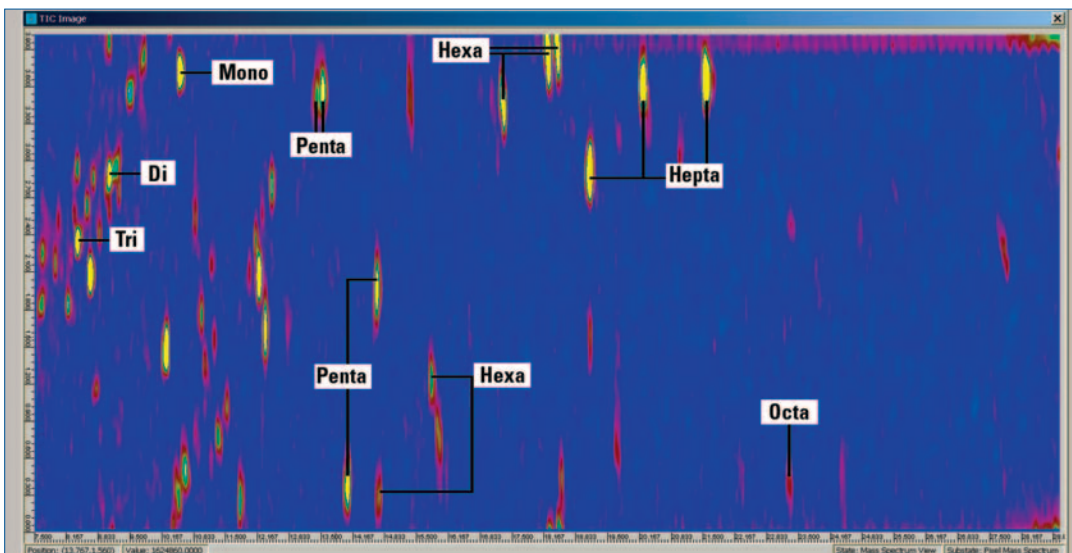
# Software

GC x GC(qMS) image of a bovine fat control sample in NCI mode

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The contour plot above shows the analysis of PCBs in bovine fat. A quadrupole mass spectrometer was used as detector after the gas chromatographic separation. In this complex sample excellent selectivity for the PCBs was achieved using the quadrupole MS in NCI mode (Negative Chemical Ionization).

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**Shimadzu Europa GmbH**